


# Badger Docs - Enable and Disable

Before a device may be used you must Enable it in Badger. Most equipment has a physical interlock which will prevent some key component from functioning unless the device is enabled in Badger. Unless instructed to do so by staff a member should never enable equipment for use by other members! Members must be “qualified” by lab staff prior to having the ability to enable a device. When you have been qualified by staff it may take a few minutes before the qualification propagates out to the system. You may need to log off and back on to get any new qualifications. Charges accrue from the moment a device is enabled and continue until the device is disabled.

## Enable a Device –

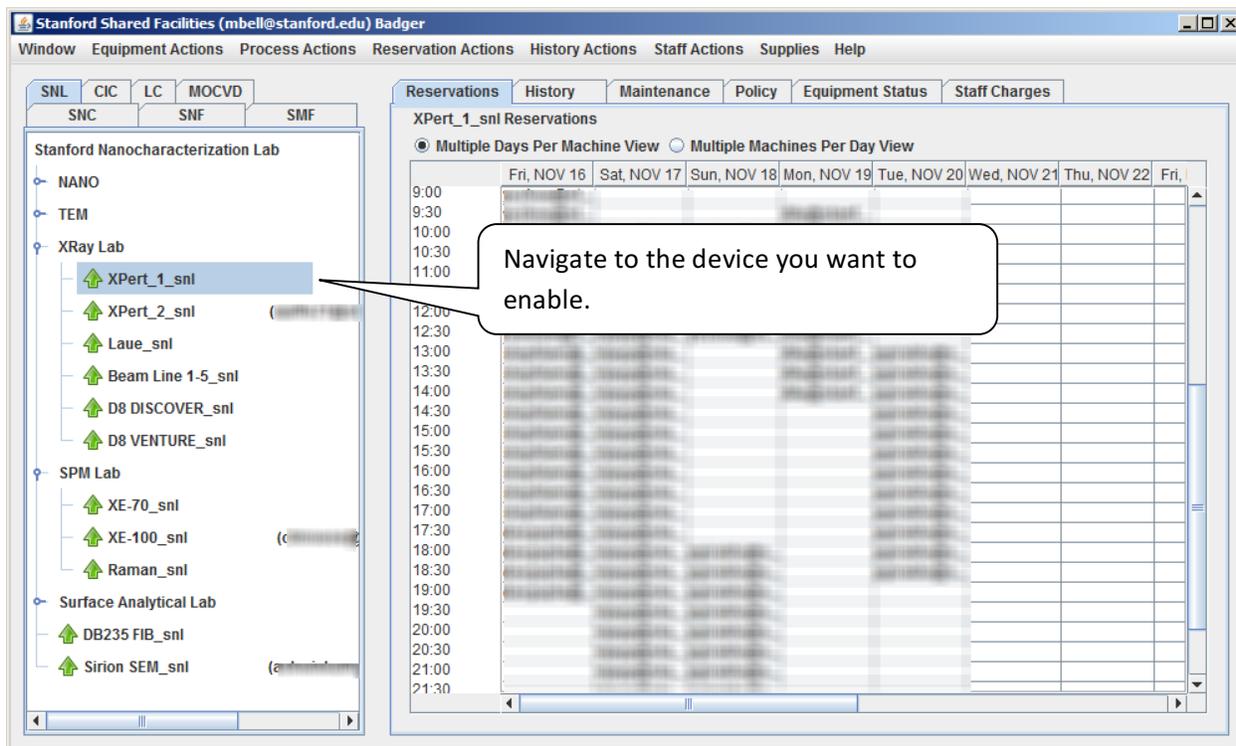

# Badger Docs - Enable and Disable

Stanford Shared Facilities (mbell@stanford.edu) Badger

Window Equipment Actions Process Actions Reservation Actions History Actions Staff Actions Supplies Help

SNL  
SN  
Stanford  
NAN  
TEM  
XRd  
XPERT\_1\_snl  
XPERT\_2\_snl  
Laue\_snl  
Beam Line 1-5\_snl  
D8 DISCOVER\_snl  
D8 VENTURE\_snl  
SPM Lab  
XE-70\_snl  
XE-100\_snl  
Raman\_snl  
Surface Analytical Lab  
DB235 FIB\_snl  
Sirion SEM\_snl

Reservations History Maintenance Policy Equipment Status Staff Charges

XPERT\_1\_snl Reservations

Multiple Days Per Machine View Multiple Machines Per Day View

	NOV 18	Mon, NOV 19	Tue, NOV 20	Wed, NOV 21	Thu, NOV 22	Fri,
12:00						
12:30						
13:00						
13:30						
14:00						
14:30						
15:00						
15:30						
16:00						
16:30						
17:00						
17:30						
18:00						
19:00						
19:30						
20:00						
20:30						
21:00						
21:30						

Select Enable under Equipment Actions.

When a device is enabled you will see a member ID next to the device. You can drag the bar at the bottom of the screen to see the entire ID.

Machine Enable Information

Enabling equipment for other members may result in expulsion from the lab.

Agent support@badgerlm

Machine FEI Nova-Nano SEM

Member support@badgerlm

Project

Account


Process default

OK Cancel

The Machine Enable Information screen gives you the ability to select project, account and process information. In most cases this doesn't change.

# Badger Docs - Enable and Disable

When you are done using the machine you must disable it. If you do not disable the tool, you will continue to be charged.

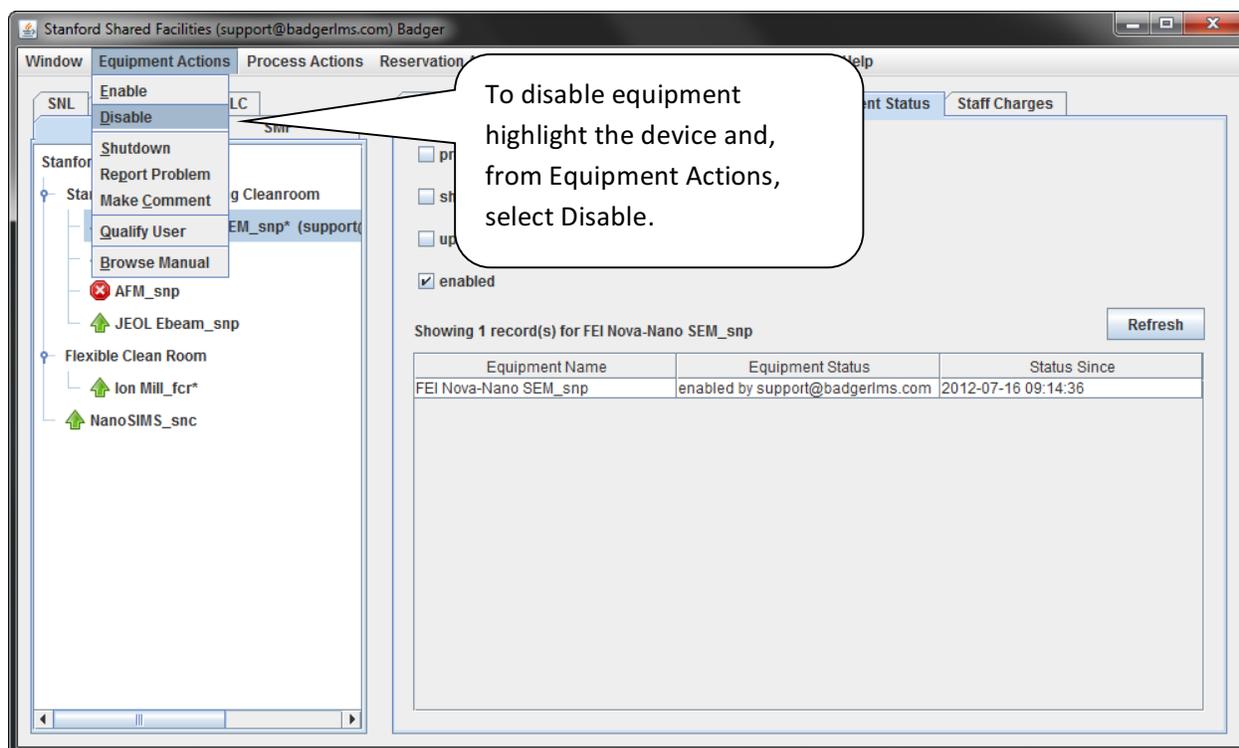

The screenshot shows the 'Stanford Shared Facilities (support@badgerlms.com) Badger' application window. On the left, a tree view shows a hierarchy of equipment categories: SNL, Stanford, and Flexible Clean Room. Under 'Stanford', the 'AFM\_snp' item is highlighted, and the 'Equipment Actions' menu is open, with 'Disable' selected. A callout box points to this menu with the text: 'To disable equipment highlight the device and, from Equipment Actions, select Disable.' The main panel shows a table with one record for 'FEI Nova-Nano SEM\_snp' with status 'enabled by support@badgerlms.com' and a 'Refresh' button.

Equipment Name	Equipment Status	Status Since
FEI Nova-Nano SEM_snp	enabled by support@badgerlms.com	2012-07-16 09:14:36

# Badger Docs - Enable and Disable

