


Badger Docs - Enable and Disable

Before a device may be used you must Enable it in Badger. Most equipment has a physical interlock which will prevent some key component from functioning unless the device is enabled in Badger. Unless instructed to do so by staff a member should never enable equipment for use by other members! Members must be “qualified” by lab staff prior to having the ability to enable a device. When you have been qualified by staff it may take a few minutes before the qualification propagates out to the system. You may need to log off and back on to get any new qualifications. Charges accrue from the moment a device is enabled and continue until the device is disabled.

Enable a Device –

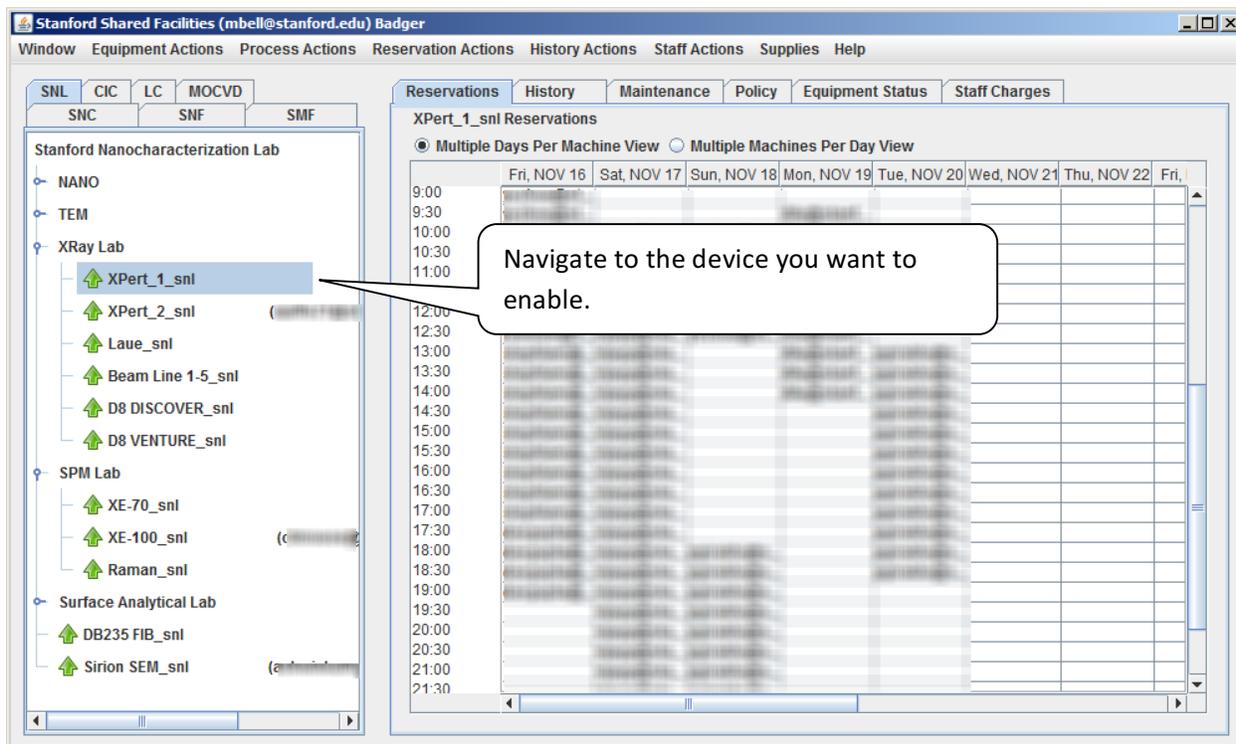


Badger Docs - Enable and Disable

The screenshot shows the 'Stanford Shared Facilities (mbell@stanford.edu) Badger' application window. The 'Equipment Actions' menu is open, with 'Enable' selected. A callout box points to this menu item with the text: "Select Enable under Equipment Actions." The main window displays a reservation table for 'XPert_1_snl' with columns for dates from NOV 18 to NOV 22 and times from 12:00 to 21:30. A second callout box points to the bottom of the reservation table with the text: "When a device is enabled you will see a member ID next to the device. You can drag the bar at the bottom of the screen to see the entire ID." The left sidebar shows a tree view of equipment categories including SNL, XRD, SPM Lab, and Surface Analytical Lab.

The 'Machine Enable Information' dialog box is shown. It contains the following fields: Agent (support@badgerlm), Machine (FEI Nova-Nano SEM), Member (support@badgerlm), Project (dropdown menu), Account (dropdown menu), and Process (default). There are 'OK' and 'Cancel' buttons at the bottom. A callout box points to the dialog with the text: "The Machine Enable Information screen gives you the ability to select project, account and process information. In most cases this doesn't change."

Badger Docs - Enable and Disable

When you are done using the machine you must disable it. If you do not disable the tool, you will continue to be charged.

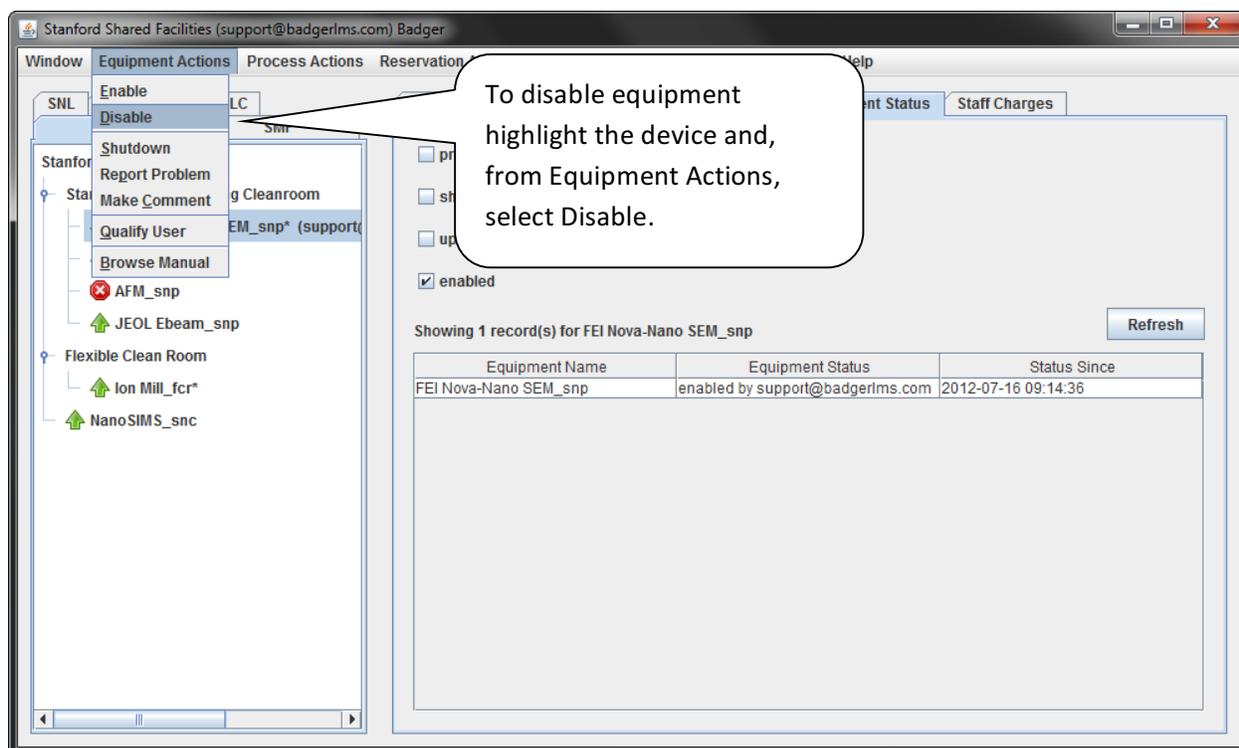

The screenshot shows the 'Stanford Shared Facilities (support@badgerlms.com) Badger' application window. On the left, a tree view shows the hierarchy of equipment, with 'AFM_snp' selected. The 'Equipment Actions' menu is open, and 'Disable' is highlighted. A callout box points to the 'Disable' option with the text: 'To disable equipment highlight the device and, from Equipment Actions, select Disable.' The main panel shows a table with one record for 'FEI Nova-Nano SEM_snp'.

Equipment Name	Equipment Status	Status Since
FEI Nova-Nano SEM_snp	enabled by support@badgerlms.com	2012-07-16 09:14:36

Badger Docs - Enable and Disable

